



Empowering the Efficient Discovery of Ultralong CDRH3 Antibodies with High-Throughput xPloration® Workflows

John J. Kim, Jonathan Theile, Katrina Yoger, Karen Padilla, Matthew Lee, Ingrid Attinost, Robert Chowdhury, David Mowrey, Soheil Khan, Lukas Fluitt, Kathryn H. Ching, Julia Noodleman, Christine N. Vuong, William D. Harriman, Philip A. Leighton, Bob Chen

OmniAb, Inc., Emeryville, CA

Abstract

The discovery of ultralong CDRH3 antibodies presents a promising frontier in therapeutic development, offering access to cryptic epitopes beyond the reach of conventional antibodies. OmniUltra™, a transgenic chicken platform engineered to express bovine-like ultralong CDRH3 antibodies in a human scaffold, combines the evolutionary advantages of avian immunology with the structural uniqueness of the stalk-knob architecture. To accelerate discovery, OmniAb® leverages xPloration®, an AI-powered, high-throughput single B-cell screening platform capable of analyzing millions of antibody-secreting cells and recovering thousands of unique sequences within hours. xPloration features a fluidics-free microcapillary system and precision laser recovery, enabling rapid capture of rare, high-value antibodies.

xPloration®: Simplifying Antibody Screening



xPloration®



How it works

1 Load

Millions of B-cells and assay reagents such as target cells or beads are easily and quickly loaded onto xPloration's microcapillary array in less than 15 min.

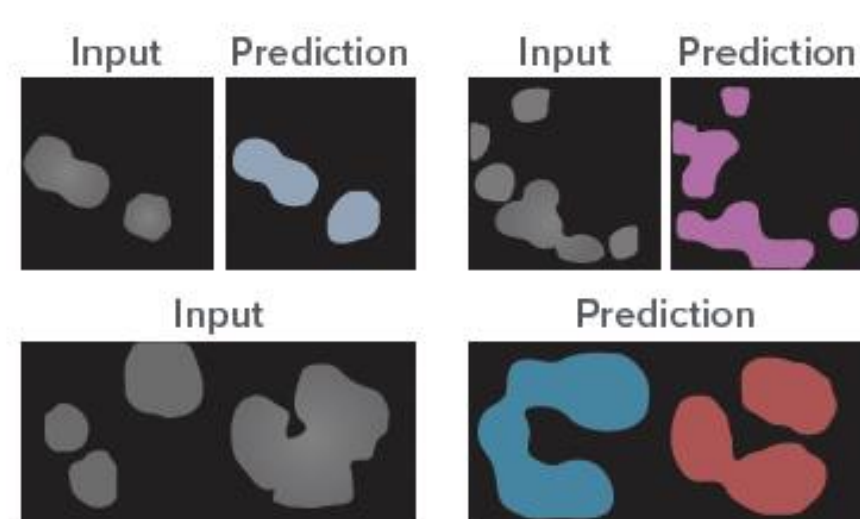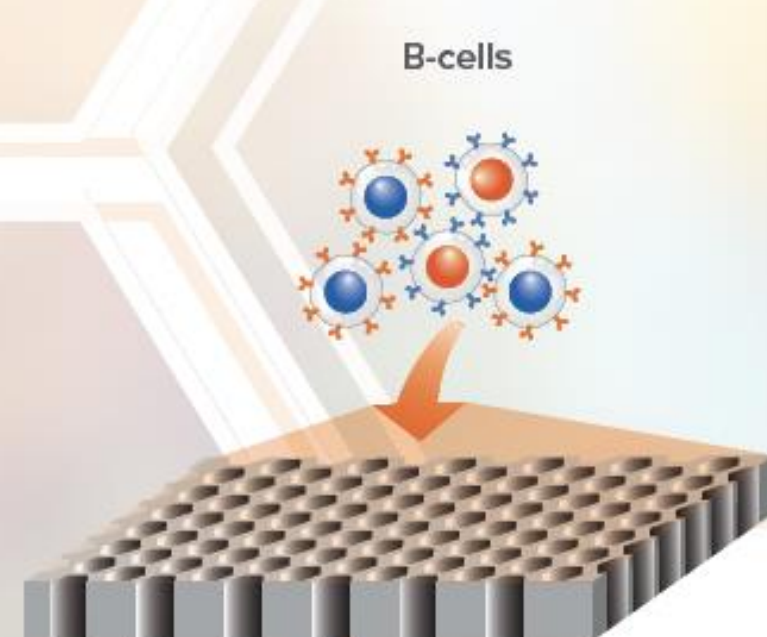


2 Analyze

xPloration instrument scans the array using AI-based image analysis to analyze every cell and identify thousands of positive hits based on your desired binding profile in real time.


3 Recover

A proprietary, precision laser recovery method gently and rapidly extracts your chosen live cells into a proprietary 96-well recovery plate, ready for immediate sequencing.

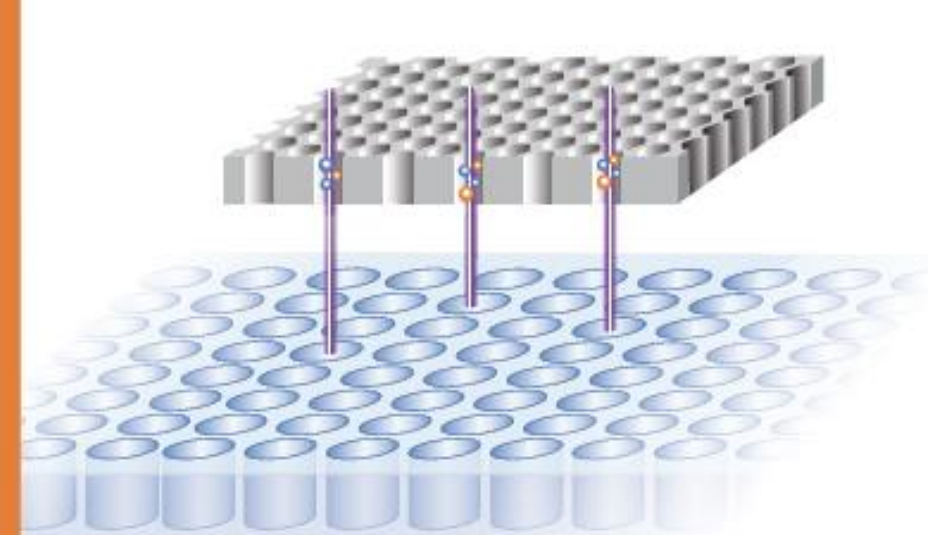


High-throughput Discovery Workflow

Case Study: PSMA is a protein highly expressed on prostate cancer cells, making it a key target for diagnosis and advanced treatments such as radioligand therapy and antibody-drug conjugates.

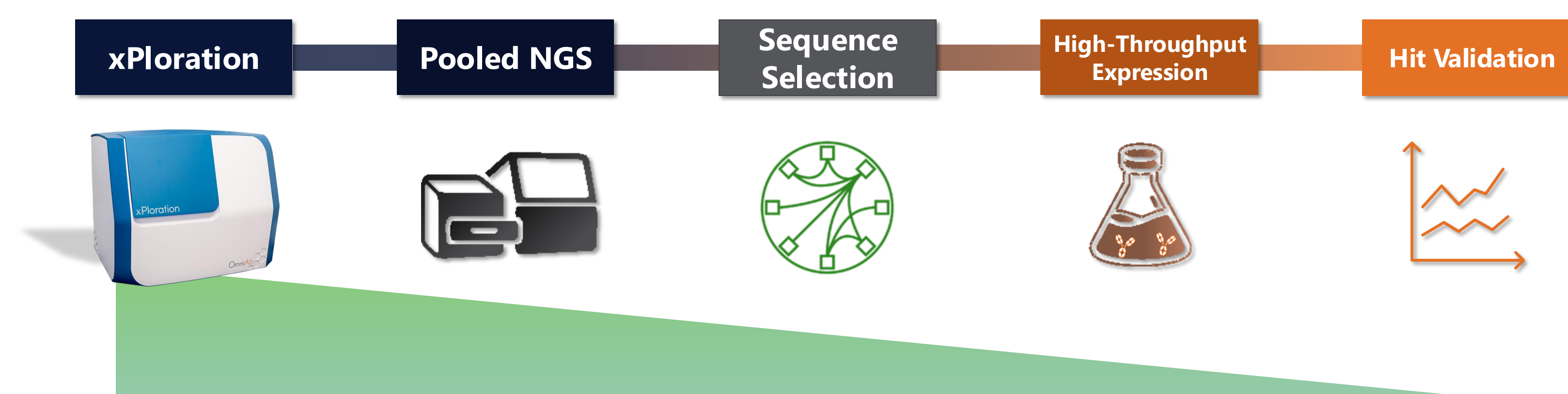

xPloration Screening

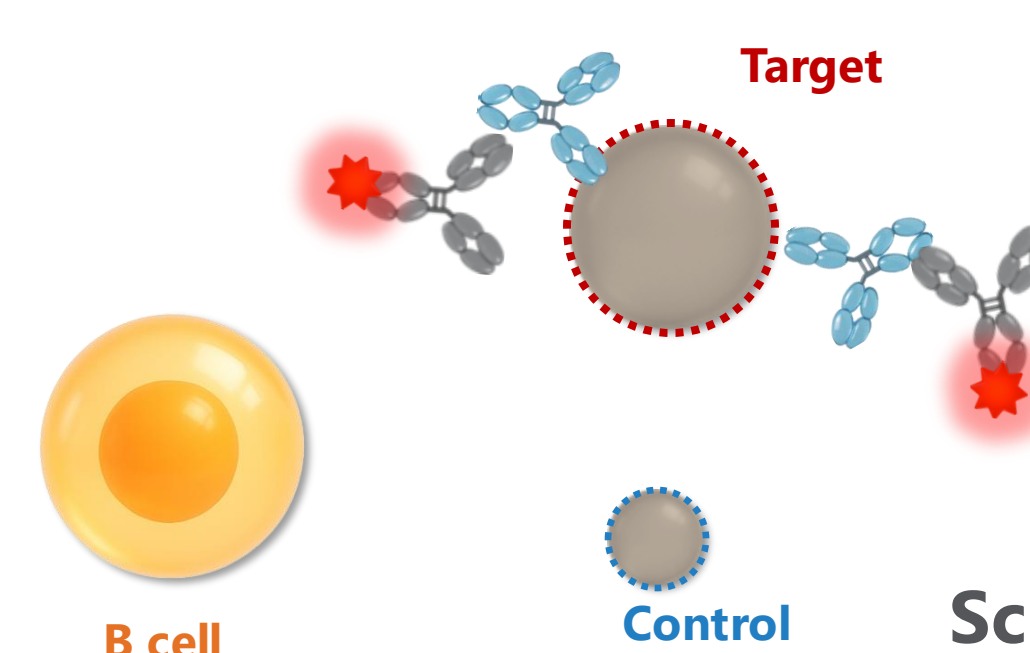


Bird	# Cells Screened	# Hits
1	14.4 M	790
2	7.2 M	1000
3	4.8 M	916
Total	26.4 M	2,706

Screening Method:

- Each bird was screened with a multiplex, protein-based, bead assay
- Hits were recovered in a pooled manner, recovering hits into 1 well




Sequence Selection

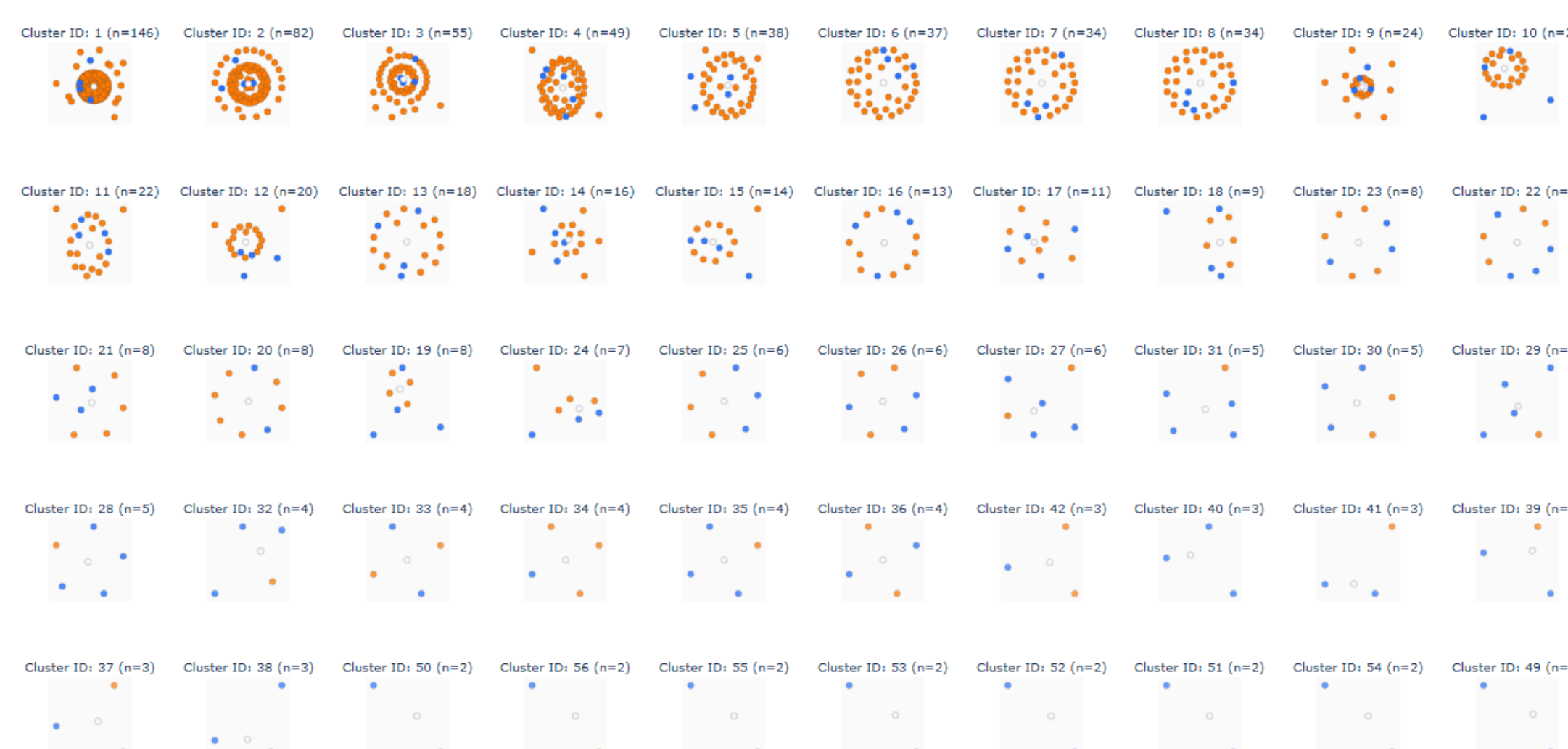

Ultralong clustering method:

- All sequences will have
 - Same Cysteine pattern
 - ≤ 3aa edit distance in knob (complete linkage)
- Selected 196 sequences
- Covered 81 clusters

Hit Validation

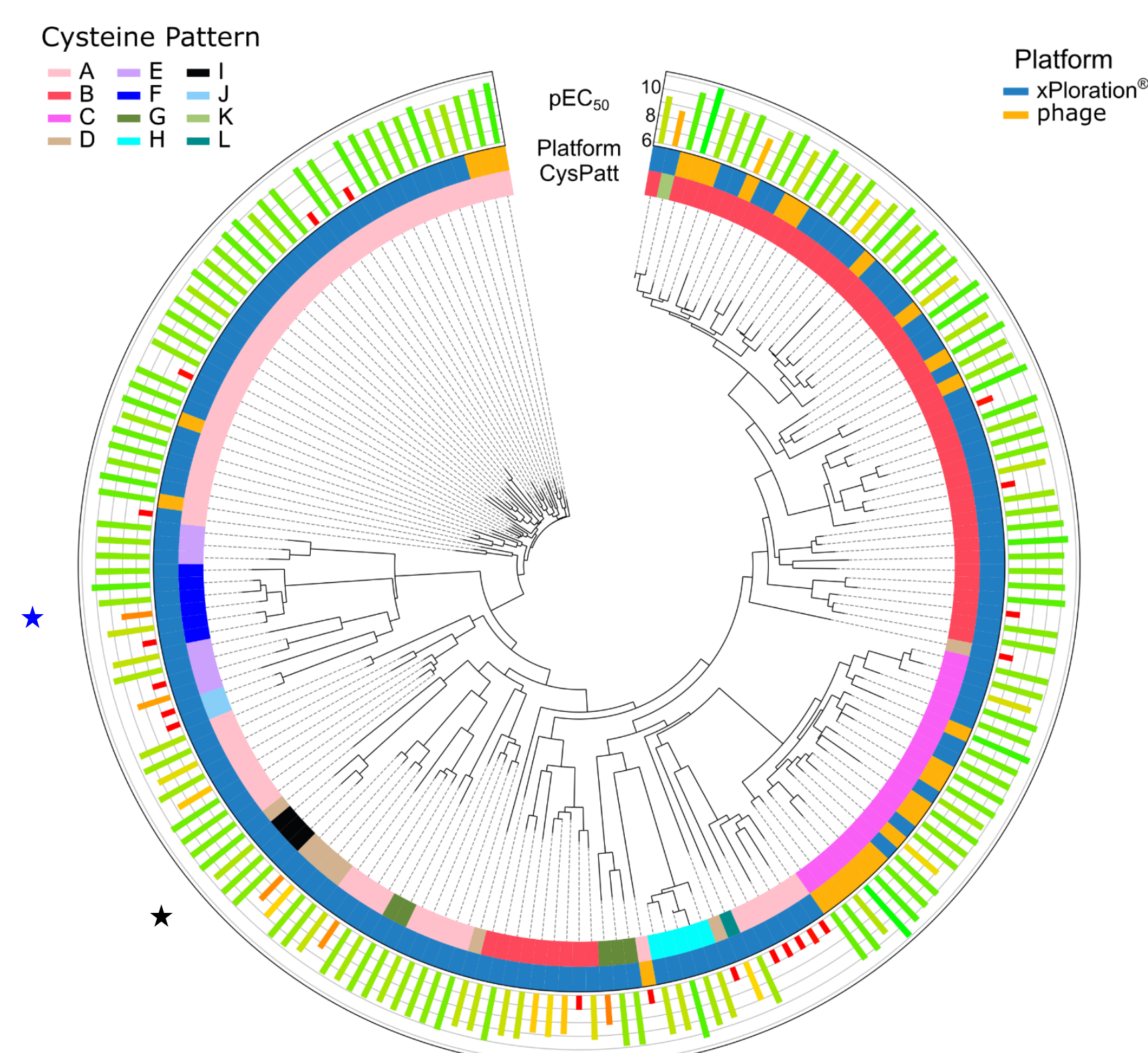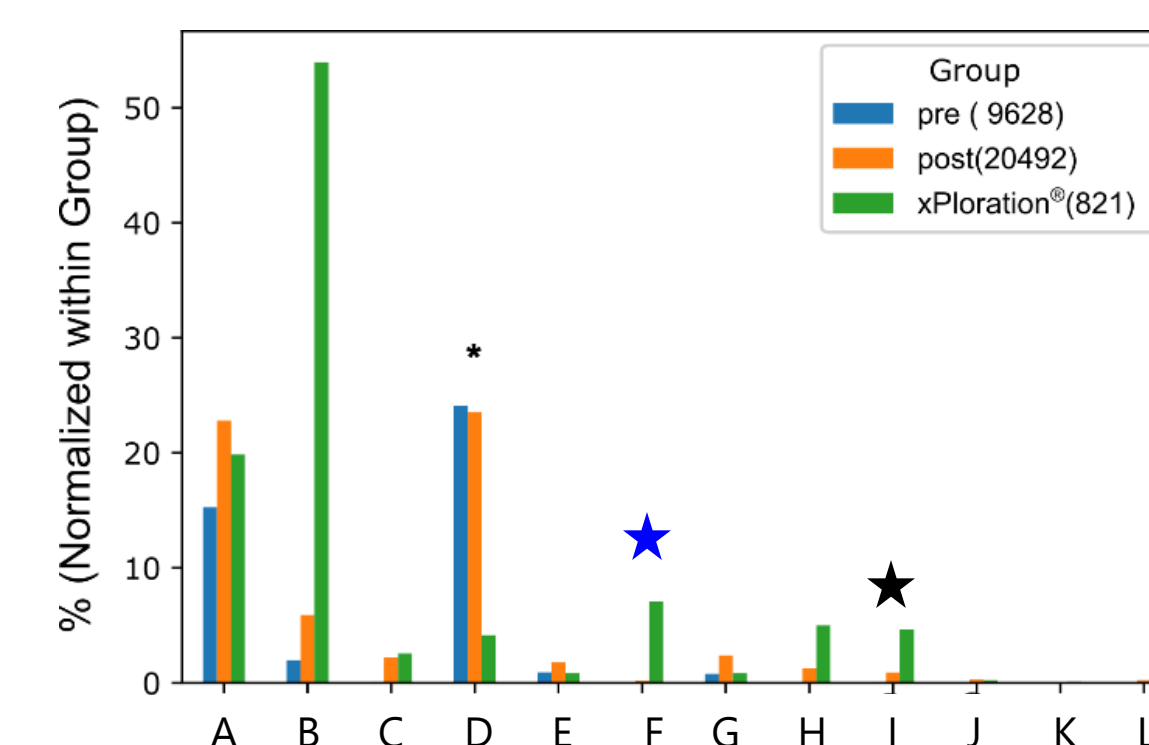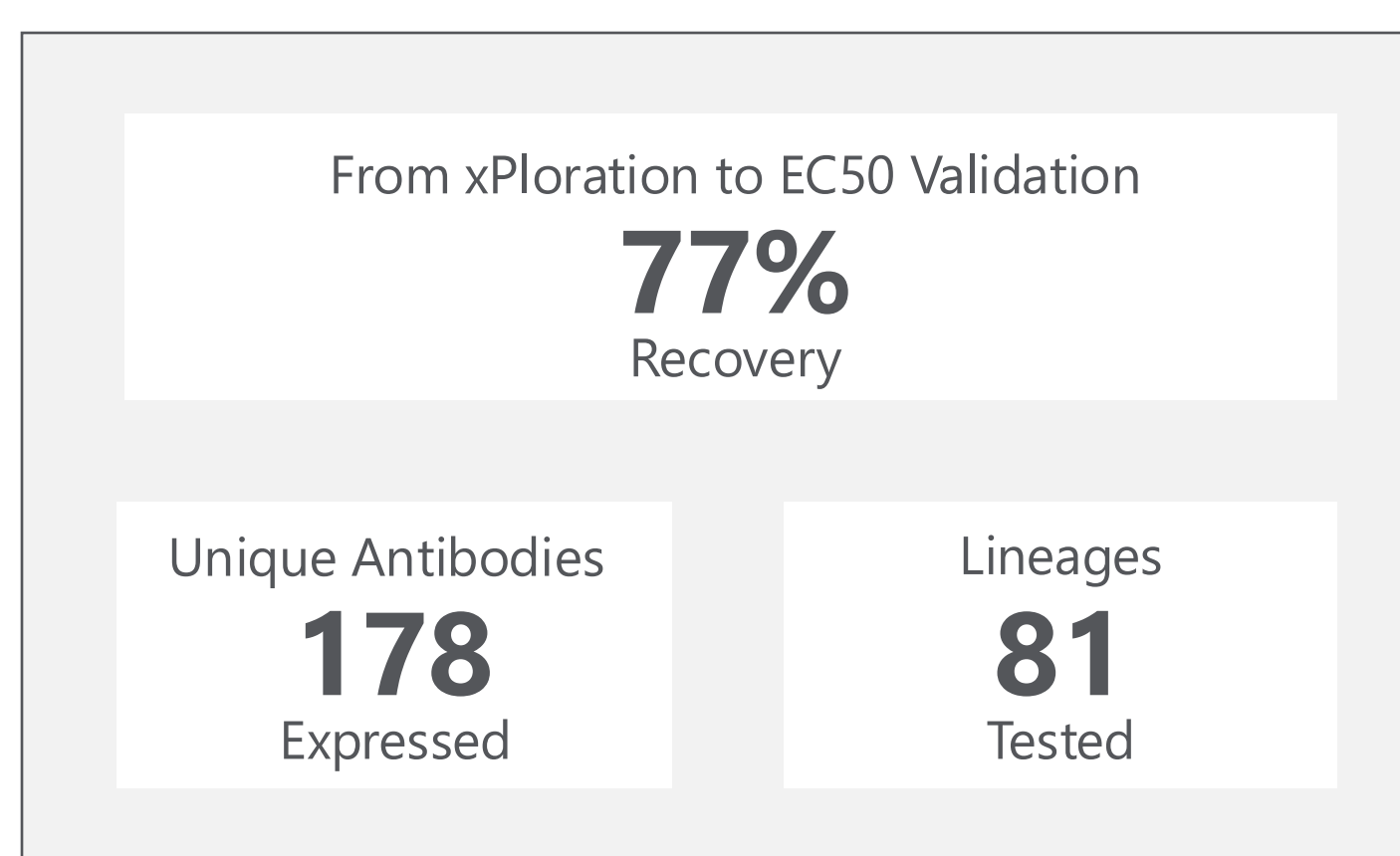

- Large range of hits from xPloration exceed phage panning
- OmniUltra™ can generate cysteine patterns distinct from germline with relevant activities against target

Conclusions

- Screening millions of cells yielded a pooled set of thousands of hits. From this repertoire, we tested and validated hundreds of candidates across more than 80 lineages.
- The combined power of xPloration® high-throughput screening and sorting, NGS, and OmniUltra™ enables large-scale discovery of antibodies with ultralong CDRH3s—and the scaffolded peptides they support.